

IL-11: IL-11R [Biotinylated] Inhibitor Screening ELISA Kit

Pack Size: 96 tests

Catalog Number: EP-159

IMPORTANT: Please carefully read this manual before performing your experiment.

For Research Use Only. Not For Use In Diagnostic Or Therapeutic Procedures

HTTP://WWW.ACROBIOSYSTEMS.COM



INTENDED USE

This kit is designed for screening of inhibitors of binding between human IL-11 and human IL-11R. It is intended for research use only (RUO).

PRINCIPLE OF THE ASSAY

This inhibitor screening ELISA kit is designed to facilitate the identification and characterization of new IL-11

pathway inhibitors. The assay takes advantage of our in house-developed binding of biotinylated human IL-11R to

immobilized human IL-11 in a functional ELISA assay, and employs a simple colorimetric ELISA platform. Briefly,

we provide you with a human Biotinylated IL-11R protein, a human IL-11 protein, an anti-IL-11 neutralizing antibody

(as method verified Std.), and Streptavidin-HRP reagent.

Your experiment will include 4 simple steps:

- 1) Coat the plate with human IL-11.
- 2) Add your molecule of interest to the tests.

3) Add human IL-11R-Biotin to bind the coated human IL-11.

4) Add Streptavidin-HRP followed by TMB or other colorimetric HRP substrate.

Finally, the half maximal inhibitory concentration (IC50) of your compound to IL-11: IL-11R binding will be

determined by comparing OD readings among different experimental groups.

MATERIALS PROVIDED

TABLE 1. MATERIALS PROVIDED (pls modify according to COA)



| Catalog | Components | Size (96 tests) | Format | Storage | | |
|-----------|----------------------------------|--------------------|--------|--------------------|-----------------------|--|
| EP159-C01 | High-bind Plate | 1 plate | Solid | 2-8°C | | |
| EP159-C02 | Human IL-11 | 55 µg | Powder | 2-8°C | -70°C after | |
| EP159-C03 | Biotinylated Human IL-11R | 20 µg | Powder | 2-8°C | reconstitution, avoid | |
| EP159-C04 | Anti-IL-11 Neutralizing Antibody | 20 µg | Powder | 2-8°C | freeze-thaw cycles | |
| EP159-C05 | Streptavidin-HRP | 5 µg | Powder | 2-8°C, avoid light | neeze thaw eyeles | |
| EP159-C06 | Coating Buffer | 12 mL | Liquid | 2-8°C | | |
| EP159-C07 | 10xWashing Buffer | 50 mL | Liquid | 2-8°C | | |
| EP159-C08 | Blocking Buffer | 50 mL | Liquid | 2-8°C | | |
| EP159-C09 | Substrate Solution | 12 mL | Liquid | 2-8°C, avoid light | | |
| EP159-C10 | Stop Solution | 7 mL | Liquid | 2-8°C | | |

REAGENTS/EQUIPMENT NEEDED BUT NOT SUPPLIED

Single or dual wavelength microplate reader with 450 nm/630nm filter ;

Centrifuge;

37 °C Incubator;

Single channel or multichannel pipettes with 10 $\mu L, 200 \; \mu L$ and 1000 μL precision ;

10 $\mu L,$ 200 μL and 1000 μL pipette tips ;

Test Tubes;

Graduated cylinder;

Deionized or distilled water for dilution ;

STORAGE AND VALIDITY INSTRUCTIONS

Unopened kit should be stored at 2°C -8°C upon receiving. Find the expiration date on the outside packaging and do not use reagents past their expiration date.

The kit should be stored as TABLE 1 after the reconstitution of lyophilized materials. The shelf life is 30 days from the date of opening.



Note:

a. Do not use reagents past their expiration date.

b. Find the expiration date on the outside packaging.

REAGENT PREPARATION

1. Restore all reagents and samples to room temperature (20-25°C) before use.

2. Reconstitute the provided lyophilized materials to stock solutions with sterile deionized water as recommended in

Tab.2, Solubilize for 15 to 30 minutes at room temperature with occasional gentle mixing. Avoid vigorous shaking or

vortex. The reconstituted stock solutions should be stored at -70°C. Avoid freeze-thaw cycles.

Note: Streptavidin-HRP stock solution should be protected from light.

TABLE 2. RECONSTITUTION METHODS FOR 96 TESTS

| Catalog | Components | Amount | Stock Solution Con. | Reconstitution Buffer and Vol. |
|-----------|----------------------------------|--------|---------------------|--------------------------------|
| EP159-C02 | Human IL-11 | 55 µg | 550 μg/mL | 100 µL, water |
| EP159-C03 | Biotinylated Human IL-11R | 20 µg | 100 μg/mL | 200 µL, water |
| EP159-C04 | Anti-IL-11 Neutralizing Antibody | 20 µg | 100 μg/mL | 200 μL, water |
| EP159-C05 | Streptavidin-HRP | 5 µg | 50 µg/mL | 100 μL, water |

RECOMMENDED PROTOCOL

1. Working solution preparation

1.1 Preparation of 1×Washing Buffer:

Dilute 50 mL 10×Washing Buffer with ultrapure water/deionized water to 500 mL.

1.2 Preparation of Dilution Buffer:

10 mL Blocking Buffer (EP159-C08) add 30 mL 1×Washing Buffer.

2. Coating

1)Dilute Human IL-11 stock solution (550 μ g/mL) to 5 μ g/mL with Coating Buffer to make Human IL-11 working

solution.

2)Add 100 μ L of Human IL-11 working solution (5 μ g/mL) to each well and leave a couple of wells uncoated for

Web: http://www.acrobiosystems.com



No-Coating Control, seal the plate with microplate sealing film and incubate overnight (or 16 hours) at 4°C.

3. Washing

Remove the remaining solution by aspiration, add 300 μ L of 1×Washing Buffer to each well, gently tap the plate for 1 minute, remove any remaining 1×Washing Buffer by aspirating or decanting, invert the plate and blot it against paper towels. Repeat the washing step above for three times.

Note: For best results, the complete removal of the Human IL-11 solution is essential. The use of a manifold dispenser or an autowasher may be necessary.

4. Blocking

Add 300 µL Blocking Buffer to each well, seal the plate with microplate sealing film and incubate at 37 °C for 1.5 hours.

5. Washing

Repeat step 3. At the same time, you can start to prepare your samples.

6. Add Samples

1)Make serial dilution of the samples as appropriate.

2)If you intend to use the provided Anti-IL-11 Neutralizing Antibody as a reference (Std.), you may dilute the antibody as recommended in Figure 1.

3)Add 50 µL of sample solution to each well according to our recommendation (Figure 2) or your own plate setup.

4)For No-Coating Control wells, please add 50 µL Dilution Buffer.

7.Binding

1) Dilute Biotinylated Human IL-11R stock solution (100 µg/mL) to 0.1 µg/mL with Dilution Buffer to make

Biotinylated Human IL-11R working solution.

2) For No-binding control wells, please add 50 μL Dilution Buffer.

3) For all other wells, please add 50 μL Biotinylated Human IL-11R working solution to the wells and mix the

samples by gently tapping the plate. Seal the plate with microplate sealing film and incubate at 37°C for 1 hour.

Note: The working solution should be prepared immediately before use and should not be stored.

FIG.1 PREPARATION OF 1:1 SERIAL DILUTIONS OF THE Anti-IL-11 Neutralizing Antibody




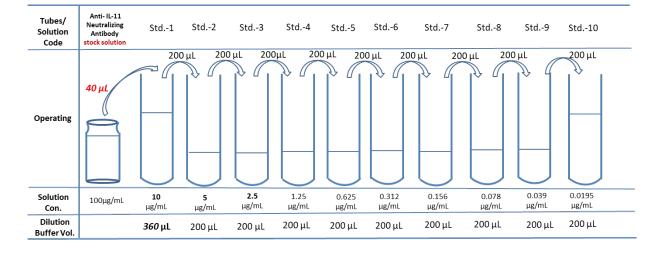


FIG.2 PLATE LAYOUT

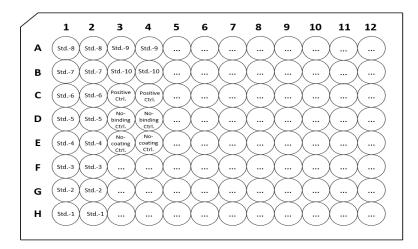

8.Washing

Repeat step 3.

9.Add Streptavidin-HRP

US and Canada:





1)Dilute Streptavidin-HRP stock solution (50 µg/mL) to 0.1 µg/mL with Dilution Buffer to make Streptavidin-HRP working solution.

2)For all wells, add 100 µL Streptavidin-HRP working solution, seal the plate with microplate sealing film and incubate at 37°C for 1 hour, avoid light.

10.Washing

Repeat step 3.

11.Substrate Reaction

Add 100 µL Substrate Solution to each well. Seal the plate with microplate sealing film and incubate at 37°C for

20 minutes. Avoid light.

12.Termination

Add 50 µL Stop Solution to each well, and gently shake the plate to allow thorough mixing.

Note: the color in the wells should change from blue to yellow.

13.Data Recording

Read the absorbance at 450 nm using UV/Vis microplate spectrophotometer.

Note: Subtracting the value read at OD_{450nm} with OD_{630nm} can be used to reduce the background noise.

SIMPLIFIED PROTOCOL

TABLE. 3 ASSAY PROTOCOL

| Steps Code | Steps | Reagents & Instruments | Reaction Conditions | Samples | No-binding Ctrl. | No-coating Ctrl. | Positive Ctrl. |
|------------|------------------------------|---------------------------------|---------------------|---------|---------------------|---------------------|-------------------|
| 1 | Working fluid preparation | N/A | N/A | N/A | N/A | N/A | N/A |
| 2 | Coating | Human IL-11 Working Solution | 4°C for overnight | 100 µL | 100 µL | | 100 µL |
| 3 | Washing | 1XWash Buffer | Wash for 3 times | 300 µL | 300 µL | 300 µL | 300 µL |
| 4 | Blocking | Blocking Buffer | 37°C for 1.5 hours | 300 µL | 300 µL | 300 µL | 300 µL |
| 5 | Washing | 1XWash Buffer | Wash for 3 times | 300 µL | 300 µL | 300 µL | 300 µL |
| 6 | Add Samples | Samples | | 50 µL | | _ | _ |
| | | Dilution Buffer | _ | _ | 50 µL | 50 µL | 50 µL |

US and Canada:

Tel: +1 800-810-0816

Web: http://www.acrobiosystems.com

7 /9

Asia and Pacific:

Tel: +86 400-682-2521

E-mail: order@acrobiosystems.com



| 7 | Binding | Biotinylated Human IL-11R Working Solution | Mix by gentle tapping, incubate at 37°C for 1 hours | 50 µL | | 50 µL | 50 µL |
|----|--------------------|---|--|--------|--------|--------|--------|
| | | Dilution Buffer | | _ | 50 µL | | — |
| 8 | Washing | 1XWash Buffer | Wash for 3 times | 300 µL | 300 µL | 300 µL | 300 µL |
| 9 | Streptavidin-HRP | Streptavidin-HRP Working Solution | 37°C for 1 hours | 100 µL | 100 μL | 100 µL | 100 µL |
| 10 | Washing | 1XWash Buffer | Wash for 3 times | 300 µL | 300 µL | 300 µL | 300 µL |
| 11 | Substrate Reaction | Substrate Solution | 37°C for 20 minutes | 100 µL | 100 µL | 100 µL | 100 µL |
| 12 | Termination | Stop Solution | Mix by gentle tapping | 50 µL | 50 µL | 50 µL | 50 µL |
| 13 | Data Recording | UV/Vis spectrophotometer | Measure absorbance at 450 nm, with the correction wavelength set at 630 nm | | | | |

Note for TAB. 3:

- 1) Samples: Your samples of interest.
- 2) No-binding Ctrl.: Reaction without Biotinylated Human IL-11R added. The absorbance should be around 0.05(< 0.1) at 450 nm.
- 3) No-coating Ctrl.: Reaction without Human IL-11 coated on the wells. The absorbance should be around 0.05 (< 0.1) at 450 nm.
- 4) Positive Ctrl.: Determined the max value in 450nm absorbance, when out of inhibitors.
- 5) It is recommended that all samples, controls and standards should be done in duplicates.

PRECAUSIONS

- 1. This kit is for research use only and is not for use in diagnostic or therapeutic applications.
- 2. This kit should be used according to the provided instructions.
- 3. Do not mix reagents from different lots.
- 4. All reagents should be balanced to room temperature (20°C-25°C) before use.
- 5. This kit should be stored at 2°C-8°C.
- 6. Please prepare the working solution of each component according to the needs of the experiment. Except for

1x Washing Buffer, all prepared working solution is for one-time use and cannot be stored.


METHOD VERIFICATION

INHIBITION OF IL-11: IL-11R [Biotinylated] BINDING BY ANTI-IL-11 NEUTRALIZING ANTIBODY

Web: http://www.acrobiosystems.com



Serial dilutions of Anti-IL-11 Neutralizing antibody (Catalog # EP159-C04) (1:1 serial dilution, from 10 µg/mL to 0.0195µg/mL) was added into IL-11: Biotinylated IL-11R binding reactions. The assay was performed according to the protocol described below. Background was subtracted from data points prior to log transformation and curve fitting.

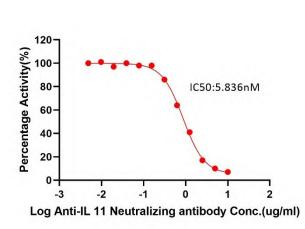

| Anti-IL-11R Neutralizing Antibody(µg/ml) | Anti- IL-11R Neutralizing Antibody.(nM) | Mean Abs.(OD450) | Percentage Activity(%) |
|--|---|---------------------|---------------------------|
| 0.000 | 0.000 | 2.685 | 100% |
| 0.005 | 0.033 | 2.694 | 100% |
| 0.010 | 0.065 | 2.700 | 100% |
| 0.020 | 0.130 | 2.612 | 97% |
| 0.039 | 0.260 | 2.675 | 99% |
| 0.078 | 0.521 | 2.643 | 98% |
| 0.156 | 1.042 | 2.632 | 98% |
| 0.313 | 2.083 | 2.304 | 86% |
| 0.625 | 4.167 | 1.715 | 64% |
| 1.250 | 8.333 | 1.100 | 41% |
| 2.5 | 16.667 | 0.468 | 17% |
| 5 | 33.333 | 0.264 | 10% |
| 10 | 66.667 | 0.189 | 7% |
| No Coating | | 0.048 | |
| No Binding | | 0.052 | |

For each experiment, a standard curve needs to be set for each micro-plate, and the specific OD value may vary depending on different laboratories, testers, or equipments. The example data is for reference only